



Saturn-X Commutator

Powered Commutator for maximum subject freedom of movement

© Neuralynx, Inc. 105 Commercial Drive, Bozeman, MT 59715 Phone 406.585.4542 • Fax 866.585.1743 www.Neuralynx.com support@Neuralynx.com

Revision 1.0 02/19/2014

Table of Contents

1	Docu	ment Overview	. 4			
2	2 Saturn-X Commutator Overview					
3	What	s Included with the Saturn-X Commutator?	. 5			
	3.1 S	aturn-X Commutator	. 5			
	3.2 S	aturn-X Power Supply	. 5			
	3.3 A	dditional Items Sold Separately	. 5			
	3.3.1	TETH-XTN-MM	. 5			
	3.3.2	Tether Interconnect	. 6			
	3.3.3	FORJ Mount	. 6			
	3.3.4	ADPT-HS-36-PSR-36	. 6			
	3.3.5	ADPT-HS-18-DUAL-PSR	. 7			
	3.3.6	ADPT-HS-27-PSR-36	. 7			
	3.3.7	Remote	. 7			
	3.4 E	lectrostatic Sensitive Equipment	. 7			
4	Quick	Start	. 8			
	4.1 S	aturn-X Commutator Setup	. 8			
	4.2 S ⁻	tart Cheetah	. 9			
	4.3 D	rive Signal into the Saturn-X Commutator	. 9			
5	Comp	outer Control	13			
	5.1 Ir	stalling the USB Drivers	13			
	5.2 S	etting Up the Terminal Window	13			
	5.3 T	erminal Window Options	15			
	5.3.1	Set Min Speed	15			
	5.3.2	Set Max Speed	15			
	5.3.3	Set Remote Speed	15			
	5.3.4	Set Acceleration Speed	15			
	5.3.5	Speed Ramp	15			
	5.3.6	Set Stop Angle	15			
	5.3.7	Set Start Angle	15			
	5.3.8	Calibrate Home Position	16			
	5.3.9	LED Speed indicator	16			
	5.3.10	Reset Factory Defaults	16			
	5.3.11	Reprint Menu	16			
6	Moun	ting	17			
7	Optic	al and Fluid Integration	18			
8	Gloss	ary	19			
9	Apper	ndix	20			
	9.1 S	aturn-X Commutator LED Indicator Descriptions	20			
	9.2 S	aturn-X Commutator Dimensions	21			

List of Figures and Tables

Figure 3-1 Saturn-X Commutator	5
Figure 3-2 Saturn-X Power Supply	5
Figure 3-3 TETH-XTN-MM	5
Figure 3-4 Tether Interconnect	6
Figure 3-5 FORJ Mount	6
Figure 3-6 ADPT-HS-36-PSR-36	6
Figure 3-7 ADPT-HS-18-DUAL-PSR	7
Figure 3-8 ADPT-HS-27-PSR-36	7
Figure 3-9 Remote	7
Figure 4-1 Hardware Connections	8
Figure 4-2 Bank 1, 2, 3 and 4 Switches Up. Reference Switch Down	9
Figure 4-3 Bank 2, 3 and 4 Switches Up. Bank 1 and Reference Switches Down	. 10
Figure 4-4 Bank 3 and 4 Switches Up. Bank 1, 2 and Reference Switches Down	. 10
Figure 4-5 Bank 4 Switches Up. Bank 1, 2, 3 and Reference Switches Down	. 11
Figure 4-6 All Switches Down (1000µV)	. 11
Figure 4-7 All Switches Down (25µV)	. 12
Figure 4-8 All Switches Up (25µV).	. 12
Figure 5-1 Driver Installation Window	. 13
Figure 5-2 Found New Hardware	. 13
Figure 5-3 COM Port Number	. 14
Figure 5-4 Terminal Window Setup	. 14
Figure 5-5 Terminal Window	. 15
Figure 7-1 FORJ Integration	. 18
Figure 9-1 LED Status Indicators	. 20
Figure 9-2 Saturn-X Commutator Dimensions – Front View	. 21
Figure 9-3 Saturn-X Commutator Dimensions – Side View	. 22

1 Document Overview

This document describes the specifications and features of the Saturn-X Commutator. It also explains how to setup your Saturn-X, test it, and use it during normal operation. There is a glossary at the end of the document.

2 Saturn-X Commutator Overview

This Saturn-X Commutator is a powered slip ring commutator that allows tethers connected to the research subject to rotate independently of the data acquisition system.

Features:

- Power Commutator
- High Sensitivity
- Up to 160 channels
- Adds Minimal Addition Noise
- User Programmable
- Optional Remote Control

3 What's Included with the Saturn-X Commutator?

There are six different channel counts available the; Saturn Mini, Saturn-1, Saturn-2, Saturn-3, Saturn-4, and Saturn-5.

3.1 Saturn-X Commutator

- 1 to 5 Channel Modules
- 32 AD Channels, 4 References, 2 Stimulus Pairs per Channel Module, except for Saturn Mini, which contains 16 AD Channels, 2 References, 2 Stimulus Pairs.
- Compatible with Neuralynx Flex Tether Adapters.
- 9.5 inches to 13.5 inches tall.
- 3.4kg to 6.5kg.

3.2 Saturn-X Power Supply

- AC/DC Power Supply.
- 24V, 1.67A.



Figure 3-1 Saturn-X Commutator



Figure 3-2 Saturn-X Power Supply

3.3 Additional Items Sold Separately

3.3.1 TETH-XTN-MM

- Extension Cable that connects the Saturn-X Commutator to the Data Acquisition System.
- Standard lengths include 1 meter, 2 meters, 5 meters, and 8 meters.



Figure 3-5 IEIH-AIN-I

Saturn-X Commutator

Revision 1.0 02/19/2014

3.3.2 Tether Interconnect

• Allows multiple TETH-XTN-MMs to connect together.



Figure 3-4 Tether Interconnect

3.3.3 FORJ Mount

- Hardware for mounting rotary joint to the top of the Saturn-X Commutator.
- 17mm and 29mm Diameter Mounts available.



Figure 3-5 FORJ Mount

3.3.4 ADPT-HS-36-PSR-36

• Allows a single HS-36 to connect to the Saturn-X Commutator.



Figure 3-6 ADPT-HS-36-PSR-36

3.3.5 ADPT-HS-18-DUAL-PSR

• Allows two HS-16/18s to connect to the Saturn-X Commutator.



Figure 3-7 ADPT-HS-18-DUAL-PSR

3.3.6 ADPT-HS-27-PSR-36

• Allows a single HS-27 or HS-16/18 to connect to the Saturn-X Commutator.



Figure 3-8 ADPT-HS-27-PSR-36

3.3.7 Remote

• Allows manual control of the Saturn-X Commutator.



Figure 3-9 Remote

3.4 Electrostatic Sensitive Equipment

All Neuralynx Equipment is Electrostatic Sensitive and should be handled with appropriate measures. Always wear a static strap and use all appropriate ESD measures when handling any electronics. Please contact Neuralynx for detailed information if you have questions.

Revision 1.0 02/19/2014

4 Quick Start

The following instructions are provided to quickly setup and test your Saturn-X Commutator setup. Note, the Saturn-X Commutator should be allowed to warm up to room temperature before performing the steps in the following section.

4.1 Saturn-X Commutator Setup

The Saturn-X Commutator connections are illustrated and described below.



Figure 4-1 Hardware Connections

Connections:

- 1. Mount the Saturn-X Commutator.
- 2. Connect the 24V Saturn-X Power Supply to the connector labeled *Power* on the Saturn-X Commutator.
- 3. Connect one end of the TETH-XNT-MM Extension Cable to Data Acquisition System.
- 4. Connect the other end of the TETH-XNT-MM Extension Cable to the connector labeled *Module 1* on the Saturn-X Commutator.
- 5. Connect the Headstage to the Flex Tether on the Saturn-X Commutator.
- 6. Connect the Headstage to the Signal Mouse.
- 7. Connect the Minirator, or other signal source, to the Signal Mouse using a BNC Cable.

Revision 1.0 02/19/2014

- 8. Turn the Bank 1, 2, 3, and 4 switches on the Signal Mouse to the Signal Position (Up). Turn the Reference switch on the Signal Mouse to the Ground Position (Down).
- Set the Minirator, or other signal source, to output a 1V_{PP} Sine Wave at 100 Hz. The Signal Mouse will reduce this signal to approximately 1mV_{PP}.
- 10. Power the Saturn-X Commutator ON.
- 11. Power the Digital Lynx SX ON.

4.2 Start Cheetah

Power On the Digital Lynx SX and wait for boot cycle to complete. On the Desktop select the *Run Cheetah* shortcut to open the Cheetah Welcome Screen. Boot Cheetah with a your preferred configuration file for your new setup. In this example 32 CSCs are used.

4.3 Drive Signal into the Saturn-X Commutator

In Cheetah, select the ACQ Button to Start Acquisition. Set the Input Range for all 32 CSCs to 1000μ V and set the Reference for all 32 CSCs to Reference 1. Observe the 32 CSCs. Each should show a reduced (approximately $1mV_{PP}$) version of the Minirator output. Refer to the figure below.

N Time	Window 1: 3000 ms < CSC.
File Do	ply Aude
C1C5	
CSCI	
CSC4	
CSC5	
CSOE	
1512	
CICE	
1.110	
CSCIE	
cscii	
CSC12	
CSC13	
CSC14	
CSCH	
CSC16	
C1017	- hannanahannanahannahannahannahannahann
csett	
- 1110	anaanaanaanaanaanaanaanaanaanaanaanaana
05020	
CSC21	
CSC22	
CSC20	
CSCM	
1.1011	hannananananananananananananananananana
C11C28	
12.7	haanaaaahaanaaahaanaaahaanaahaanaahaanaahaanaahaanahaanahaanahaanaahaanaahaanaahaanaahaah
C1(2)	hannannannannannannannannannannannannann
C5C29	
CSC20	
CSCH	
C1C12	

Figure 4-2 Bank 1, 2, 3 and 4 Switches Up. Reference Switch Down.

Switch the Bank 1 Switch to the Ground Position (Down). Observe the 32 CSCs. CSCs 1-8 should now be flatlined while CSCs 9-32 still show a reduced (roughly $1mV_{PP}$) version of the Minirator output. Refer to the figure below.

N Time W	indoe 1: 5000 ms - CSC1
File Disp	lay Audio
CSC2	
CSCI	
CSC4	
CSC5	
CSCE	
6.624	
CECE	
1111	
CSCHE	
CSCI	
CSC12	
CSC13	
CSC14	
CSCH	
CSC16	
05017	
CSCHE	
	hanna an
CSC20	
C9C21	
CSC22	
CSC20	
CSCDA	
1.12219	hannaanhannaanhannaahannaahannaanaanhannaanaa
C1028	hannannahannanahannahannahannannahannann
1.00	kananananahanananananananananananananana
C5C28	hannaanahannaanhannaahannaanhannaanhannaanhannaanhannaananhannaananhannaanhannaanhannaanhannaanhannaanha
CSC29	
CSC20	
CECH	
CSC12	hammahammahammahammahammahammahammaham

Figure 4-3 Bank 2, 3 and 4 Switches Up. Bank 1 and Reference Switches Down. Switch the Bank 2 Switch to the Ground Position (Down). Observe the 32 CSCs. CSCs 1-16 should now be flatlined while CSCs 17-32 still show a reduced (roughly 1mV_{PP}) version of the Minirator output. Refer to the figure below.

TA sume of	NOON 2. JOOD ING - COCL																		6.00	
File Displ	ay Audio				_													-		
-																				
1000																				
CSCE																				
CSCF																				
CSC4																				
CSC5																				
CSCE																				
and the second																				
CICE																				
1100																				
CSCIE																				
CSCH																				
05012																				
CSC13																				
CTICLE																				
Cacin																				
CSCIR																				
CSCIE																				
0100	00000000000	000000	anna	00000	0000	AAAA	100000	0000	000000	nana	000000	00000	00000	0000	00000	000	00000	0.0.0	0000	00000
1.11.12	~~~~~	000000	00000	00000		0000	000000		000000	0000		00000						441	Y V V	000000
CSC18	www.ww	www	www	www	nn	ww	www	www	www	vvvv	www	www	mm	\sim	www	www	www	w	MM	AMM
	ANNANA ANA				NAAA															
05030	000000000000000000000000000000000000000		AAAA	AAAAA	AAAA	AAA				10000	~~~~~	AAAAA		AAAA.		1000		A A A	AAAA	AAAAA
		000000	00000			0000	000000		000000	0000		00000	00000			0000				
CSC24	~~~~~~~~~~~	000000	~~~~	VVVVV	10000	VVVV	000000	VVVV	000000	VVVVI	100000	AAAAAA	VVVVV	10000	VVVVV	0000	~~~~~	VVV	VVVV	VVVVVV
CSC22	www.ww	www	m	www	m	ww	www	www	\sim	www	www	www	mm	M	www	www	www	\sim	~~~	mm
CSC23	~~~~~	~~~~	AAAA	~~~~	AAAA	MAAN	11111	AAAA	AAAAA	1000	ΛΛΛΛΛΛ		~~~~	AAAA	~~~~	1111	٨٨٨٨		AAAA	ΛΛΛΛΛ
	0.0.0.0.0.0.0.0	00000				0.0.01	00000											0.0.0		00000
Course	~~~~~~~~~~	~~~~~	~~~~	VVVVV	0000	0000	~~~~~		000000	0000	100000	00000	00000	10000	~~~~	VVVV	~~~~	VVV	/ / / / /	00000
110228	wwwww	www	MM	MMAA	MM	ww	\sim	$\sim \sim \sim$	www	www	www	www	www	MM	www	MAA	www	M	ANN	AMMA
C1C28	AAAAAAAAA	AAAAAA	AAAA	AAAAA	AAAA	MAAN	MAAAA	AAAA	AAAAAA	MAAAA	AAAAA	AAAAA	AAAAA	AAAA	AAAAA	AAA	AAAAA	AAA	AAAA	ΛΛΛΛΛ
	AAAAAAAAA																			
C1C28	www.www	www	www	www	mm	www	www	m	www	www	www	www	mm	MM	www	www	www	Ś	ww	www
C5C29	mmm	mm	www	mm	m	ww	www	m	www	www	mm	www	mm	M	www	www	www	w	m	mm
cscae -	000000000000000000000000000000000000000	ΛΛΛΛΛΛ	AAAA	1000	MAAA	AAA	10000	1000	ΛΛΛΛΛΛ	MAAAA	AAAAAA	AAAAA	00000	ΔΔΔΔ	00000	1000	00000	AAA	AAAA	ΛΛΛΛΛ
			00000			0000		0.000								0000				
CSCH	000000000000000000000000000000000000000	0000000	00000	VVVVV	0000	VVVV	000000	VVVV	000000	0000	100000	00000	00000	10000	00000	VVVV	00000	VVV	1000	100000
CSC12	www.ww	mm	m	m	m	ww	www	www	www	www	mm	www	mm	M	m	www	mm	Ś	ww	mm
																		_		

Figure 4-4 Bank 3 and 4 Switches Up. Bank 1, 2 and Reference Switches Down. Switch the Bank 3 Switch to the Ground Position (Down). Observe the 32 CSCs. CSCs 1-24 should now be flatlined while CSCs 25-32 still show a reduced (roughly 1mV_{PP}) version of the Minirator output. Refer to the figure below.

Revision 1.0 02/19/2014

N Time V	Window 1: 5000 ms - CSC1	
File Disp	July Audie	
C8C5		
CSCF		
CSC4		
CSC5		
CSCE		
2127		
CECE		
1111		
05010		
CSCN		
C5C12		
CSCH		
CSC14		
CSCIE		
CSCIE		
01017		
CSÓIII		
-		
05020		
05021		
05022		
C1020		
CORDA		
CITCLE		
CICA		*****
C)(2)		
CICN		
CSC/I		
CSCH		www.www.www.www.www.www.
C1012	1~~~~~~~~~~~~~~~~~~~~~~~~~~~~~~~~~~~~~~	

Figure 4-5 Bank 4 Switches Up. Bank 1, 2, 3 and Reference Switches Down. Switch the Bank 4 Switch to the Ground Position (Down). Observe the 32 CSCs. CSCs 1-32 should now be flatlined. Refer to the figure below.

11							terret in the second se
File Displa	y Aud	lio					
and the second second	-						
and the second se							
	1						
C5C2							
TOTAL OF							
1000							
CSC4							
Caca							
CSCE			 	 			
21627							
F 1 5 5							
0.000							
1.1.1.1							
15010							
CSCH							
C\$C12							
01010							
Calcia							
CSC14							
_							
CSCM							
CSCIE							
1251.47							
C S C H H							
1.11.10							
22210							
CSC20							
C9C21							
00000							
Colum							
CSC28							
CPL28							
1100000							
C1C28							
CSC28							
00000							
Courses of							
CSC28							
CICH							
CSC12							

Figure 4-6 All Switches Down (1000µV).

Set the Input Range for all 32 CSCs to 25μ V. Observe the 32 CSCs. CSCs 1-32 now show the baseline noise. Each should be less than 25μ Vpp and void of any repetitive signals. Refer to the figure below.

Revision 1.0 02/19/2014



Figure 4-7 All Switches Down (25µV).

Force the Saturn-X Commutator to spin. Observe the 32 CSCs. CSCs 1-32 should still show the baseline noise. Each should be less than 25μ Vpp and void of any repetitive signals.

Switch all the Switches to the Signal Position (Up). Observe the 32 CSCs. Each should be less than 25μ Vpp and void of any repetitive signals. Refer to the figure below.

File Display Audio								
-								
C5C5								
CSCF			and the second of the second					
CSC4								
CSCS								
CSOE					والمحمد والمحمد والمحمد والمحمد والمحافظ المحاف			
CHOT								
0100								
C 5 C 10								
05014								
CSC12								
cscu								
CSC14								
CSCH		and an a product of the standard of the standard standard standard standard standard standard standard standard						
CSC16								
05017								
CSCIE	Ne fediciel industry this parts to a loss of the liter	-						
05020								
C\$C21	ange die eine beite gester beiter							
CSC22								
05023								
65654								
0.000								
Citan								
csca								
CSC28	the second and a state of a state							
CSC20								
CSCH		and head of the second s	and the second	وجارك والماسان المرجع المرجع المرجع	Made and a second state of the second state of the	College of the design of the second		
CSC12	W. and The Links of the Links of the Links		and a second descenting of the descent		Minister description to the description of the desc			

Figure 4-8 All Switches Up (25µV).

5 Computer Control

5.1 Installing the USB Drivers

Computer control over the Saturn-X Commutator is accomplished via a Virtual COM Port and a USB to Serial UART IC in the device. The drivers for the Virtual COM Port can be found at the Neuralynx website under the Software Section. By enabling this connection, the user can access control settings for the Saturn-X Commutator. These include Speed, Acceleration, Start Angle, Stop Angle, and Home Position.

Once the driver package has been downloaded the drivers for the Virtual COM Port can be found in the Folder labeled USB Drivers. Install the drivers by running the *CDMXXXXX_Setup.exe* file. The window in the figure below will appear.



Figure 5-1 Driver Installation Window

It may take a few minutes to install the drivers. When it is finished, make sure the window states: "FTDI CDM Driver Installation process completed." Once the drivers have been installed the window will close itself.

You can now attach the Saturn-X Commutator to your computer with the USB Cable. After cycling through some Found New Hardware popups the popup shown in the following figure will appear confirming the driver was installed properly.



Figure 5-2 Found New Hardware

5.2 Setting Up the Terminal Window

PuTTY, a free Terminal Window is also included in the driver package.

Before opening the Terminal Window the COM Port Number must first be found. Do this by opening the Device Manager and expanding the Ports (COM and LPT) tab. Open the

Revision 1.0 02/19/2014

Device Manager by right-clicking on the My Computer Icon followed by selecting Properties then selecting the Hardware Tab and finally clicking Device Manager. The COM Port Number will be listed after the USB Serial Port. The COM Port Number of the figure below is **COM7**.



Figure 5-3 COM Port Number

Now that the COM Port Number has been obtained, the Terminal Window can be configured. Do this by opening the File *PuTTy.exe* found in the PuTTY Folder. The window in the following figure will appear.

🞇 PuTTY Configuration		×
Category:		
Session Logging Terminal Keyboard Bell Features Window Appearance Behaviour Translation	Basic options for your PuTTY se Specify the destination you want to conne Serial line COM7 Connection type: C Raw C Ielnet C Rlogin C SSI Load, save or delete a stored session Saved Sessions	ssion ct to Speed 115200 H
Selection Colours Connection Oata Proxy Telnet Rlogin SSH Serial	Default Settings	Load Sa <u>v</u> e Delete
<u>A</u> bout	Close window on e <u>x</u> it: C Always C Never Only on o Open	lean exit <u>C</u> ancel

Figure 5-4 Terminal Window Setup

Enter the COM Port Number found earlier into the text box below Serial Line. Select Default Settings under Saved Sessions and click Save. Click Cancel when finished. The Terminal Window has now been configured.

Now open *PuTTY.exe* in the root folder. Press "99" then Enter on the Keyboard, the Terminal Window should display the Auto Mode choices as shown in the figure below. If the Terminal Window does match the figure below the Terminal Window has been configured properly.

Revision 1.0 02/19/2014

🛃 P	uTTY				↔	_ D ×			
Neu	Neuralynx Commutator Controller								
1) 2) 3) 4) 5) 6) 7) 8) 9)	Set Min Speed Set Max Speed Set Remote Speed Set Acceleration Speed Ramp Set Stop Angle Set Start Angle Calibrate home position LED Speed indicator Enable Smart Remote		10000 25000 10000 3 1 3 176 1 0						
98) 99)	Reset Factory Defaults Reprint Menu		-						
Ent	er menu option number and	press	Enter:						

Figure 5-5 Terminal Window

5.3 Terminal Window Options

5.3.1 Set Min Speed

Sets the minimum pulse rate for the Saturn-X Commutator motor.

5.3.2 Set Max Speed

Sets the maximum pulse rate for the Saturn-X Commutator motor.

5.3.3 Set Remote Speed

Sets the pulse rate for the Saturn-X Commutator motor in Manual Mode.

5.3.4 Set Acceleration Speed

When combined with Ramp Setting, sets the speed at which the Saturn-X Commutator motor accelerates from its minimum pulse rate to its maximum pulse rate, based on the amount of tether displacement.

5.3.5 Speed Ramp

When combined with Acceleration Setting, sets the speed at which the Saturn-X Commutator motor accelerates from its minimum pulse rate to its maximum pulse rate, based on the amount of tether displacement.

5.3.6 Set Stop Angle

Sets the number of displacement degrees from the home position at which the motor will stop spinning.

5.3.7 Set Start Angle

Sets the number of displacement degrees from the home position at which the motor will start spinning.

Revision 1.0	
02/19/2014	

5.3.8 Calibrate Home Position

Calibrates the home position of the Saturn-X Commutator.

5.3.9 LED Speed indicator

Enables and Disables the Status LED on the Saturn-X Commutator.

5.3.10 Reset Factory Defaults

Resets the Saturn-X Commutator to factory defaults. The Saturn-X Commutator will most likely immediately begin spinning after selecting this option.

5.3.11 Reprint Menu

Simply reprints the Neuralynx Commutator Controller Menu.

6 Mounting

The Saturn-X Commutator is designed to be mounted above an experiment location, usually at the ceiling of the experiment room. The Saturn-X Commutator platform should be mounted as parallel to the floor as possible. Vibration in the mounting structure can cause the torque sensor to shift enough to cause motor rotation. If this happens with nothing attached to provide counter torque, the motor will spin continuously until counter torque is applied or power is removed. The chance of this unintended rotation occurring can be reduced by using the sensitivity threshold jumpers.

Due to the difficulty of locating the Cheetah system near the ceiling, the Saturn-X Commutator is always used with the Neuralynx tether extension, which also protects the delicate wiring from damage in a potentially hostile ceiling environment. The Neuralynx headstage buffers are able to drive up to 25 feet of tether extension cable with no increase in system noise. In most cases, the use of the Saturn-X Commutator and extension cabling should be transparent to the performance of the system. It is suggested, however, that the power cabling for the Saturn-X Commutator be run at least half of one meter away from the tether extension (along its entire length), in order to avoid conducting power supply noise into the electrode signals. This is not a hard requirement but is strongly recommended.

The flex circuit tether exiting the bottom of the Saturn-X Commutator is connected to the headstage tether using the appropriate adapter board. The tether and headstage assembly are usually left plugged in after mounting. The animal is disconnected and reconnected to the hanging tether and headstage each day.

7 Optical and Fluid Integration

The Saturn-X Commutator comes pre-designed for optical and fluid integration. This mount can be purchased separate and can be added at anytime after initial installation. The Saturn-X Commutator unit works independently of the FORJ or Fluid Rotary Joint. The mount does not include either fluid or optical joint. These must be purchased separately.



Figure 7-1 FORJ Integration

8 Glossary

CSC – Neuralynx acronym for Continuously Sampled Channel.

EIB – Neuralynx acronym for Electrode Interface Board.

HS – Neuralynx acronym for Headstage.

Saturn-X Commutator – Powered slip ring commutator. X denotes the number of Channel Modules (ie. Saturn-1 contains one Channel Module).

9 Appendix

9.1 Saturn-X Commutator LED Indicator Descriptions



- Saturn-X Commutator powered Off.
- Saturn-X Commutator powered On.
- Power Good.
- Saturn-X Commutator powered On.
- Power Good.
- Motor Spinning.
- Saturn-X Commutator powered On.
- Power Good.
- Device in Manual Mode
- USB Cable connected to Saturn-X Commutator.
- Serious power problem.
- Contact Neuralynx for support.
- Saturn-X Commutator powered On.
- Power Good.
- Serious Saturn-X Commutator problem.
- Contact Neuralynx for support.

Figure 9-1 LED Status Indicators



9.2 Saturn-X Commutator Dimensions

10.000" Figure 9-2 Saturn-X Commutator Dimensions – Front View



11.000" Figure 9-3 Saturn-X Commutator Dimensions – Side View